

# Agilent 7700 Series ICP-MS Techniques and Operation

Course Number R1777A

Student Lecture Manual





**Agilent Technologies**

# Agilent 7700 Series ICP-MS Techniques and Operation

Course Number R1777A

Student Lecture Manual

Agilent 7700 MassHunter Workstation  
Printed in September, 2009

## **Notice**

The information contained in this document is subject to change without notice.

Agilent Technologies makes no warranty of any kind with regard to this material, including but not limited to the implied warranties of merchantability and fitness for a particular purpose.

Agilent Technologies shall not be liable for errors contained herein or for incidental, or consequential damages in connection with the furnishing, performance, or use of this material.

No part of this document may be photocopied or reproduced, or translated to another program language without the prior written consent of Agilent Technologies, Inc.

Agilent Technologies, Inc  
3750 Brookside Pkwy  
Suite 100  
Alpharetta, GA 30022

© 2009 by Agilent Technologies, Inc.

All rights reserved

Printed in the United States of America

## Table of Contents

<b>AGILENT 7700X AND 7700S ICP-MS PRINCIPLES AND OPERATION .....</b>	<b>1</b>
In This Section .....	2
ICP-MS's Broad Reach.....	3
ICP-MS Application End-Use .....	4
ICP-MS Geographical Perspective .....	5
Reasons for Popularity .....	6
Agilent Contributions to ICP-MS .....	8
Software Contributions .....	9
Training Program .....	10
Presentation Slides .....	12
Figures of Merit .....	14
Too Hard to Understand?.....	15
<b>FUNDAMENTALS OF ICP-MS AND AGILENT 7700 DESIGN FEATURES.....</b>	<b>17</b>
In This Section .....	18
Measuring Matter.....	19
Spectrometric Techniques.....	20
The Source and Analyzer.....	21
Atomic Structure.....	22
Atomic Ions.....	24
Ionization Potential .....	26
Isotopes and Abundance .....	27
Control of Ion Motion.....	28
Mass 6 to 260.....	29
Neutral Particles.....	30
Doubly Charged Ions .....	31
Ion Energy and Spread of Ion Energy.....	32
Anatomy of a Signal .....	35
The Conventional ICP-MS Mass Spectrum.....	36
Liquid Sample Introduction .....	37
Liquid-Aerosol Sample Introduction Components.....	38
Argon Gas Lines .....	39
Nebulizer Basics .....	40

Alternative Nebulizer Choices.....	41
Nebulizer Comparison .....	42
Spray Chamber.....	43
ICP Torch.....	45
Inductively Coupled Plasma .....	46
Aerosol Particle Injection into the Plasma.....	47
Interface: Sample Entering Vacuum.....	50
Key Interface Components .....	51
Ion Lenses .....	52
Quadrupole Mass Spectrometer .....	54
QP Data Acquisition .....	55
Agilent 7700 QP Mass Analyzer .....	56
Bias Voltage Applied to QP and ORS .....	57
Multiplier (EM) Basics .....	58
Multiplier Optimization .....	59
Agilent 7700 Pumping System .....	60
Agilent 7700 ICP-MS Full System View .....	61
Utilities.....	62
<b>BACKBONE OF MASSHUNTER SOFTWARE FOR AGILENT 7700 INSTRUMENT</b>	
<b>SETUP/STARTUP .....</b>	<b>63</b>
In This Section .....	64
Common File and Folder Extensions.....	65
File Naming .....	66
Desktop and Software Icons .....	68
Configuration .....	69
ALS Positioning: ASX-500 .....	74
Configuration Setup for Agilent I-AS.....	75
ALS Positioning: I-AS.....	76
Starting the Instrument.....	77
Menu Row for Instrument Control .....	78
Meters .....	79
Peristaltic Pump .....	80
Torch Box and Sample Introduction Access .....	82
Torch Box with Quartzware Assembled.....	83

ICP Ignition to ANALYSIS State.....	84
Setup Modes.....	85
Setup Mode: Mixing Tee .....	86
HMI: High Matrix Introduction .....	87
Setup Mode: Inert Sample Introduction.....	90
Setup Mode: ISIS.....	91
<b>AGILENT 7700 TUNING.....</b>	<b>93</b>
In This Section .....	94
Tuning = Optimization.....	95
Tune Modes .....	96
Primary Tuning: No-Gas Mode .....	97
Key Prerequisite Goals for Tuning .....	98
Tuning Parameters .....	99
Tuning Solutions.....	100
Tuning Parameter Adjustment Overview .....	101
“Stability” Is a Matter of Timeframe .....	102
Main Tuning Window.....	103
Main Tune Window: Acq. Parameters.....	105
System Wide Parameters .....	106
Automation of Tuning.....	107
Autotune Software Feature .....	108
Plasma Conditions Auto Selection.....	109
Autotune: Robust Tuning.....	110
Autotune: User Defined .....	111
Autotune Activity Window.....	112
Manual Tuning (No-Gas Mode): Sensitivity .....	113
Tuning Goal: Good Sensitivity at Each Mass.....	114
Manual Tuning (No-Gas Mode): Stability.....	115
Tuning (No-Gas Mode): Robust = Low Oxides .....	116
Oxides and Nebulizer Gas Flow .....	117
Common No-Gas Mode Tune Settings.....	118
QP Tuning: Resolution/Axis Window .....	119

EM Tuning .....	121
Detector Tuning: Pulse/Analog (P/A).....	122
P/A Tune Report .....	123
Saving Tune Files .....	124
Tune Report .....	125
Review: How Optimization is Measured.....	126
Manual Optimization: Know Your Tuning Goal.....	127
Tuning Verification Summary .....	128
Significance of the Tune File.....	130
<b>ICP-MS INTERFERENCES.....</b>	<b>131</b>
In This Section .....	132
ICP-MS Interferences .....	133
What Is a Matrix?.....	134
Impact of Dissolved Solids .....	135
Suppression Effects.....	136
Space Charge Effects .....	137
ICP-MS Nebulization Effects .....	138
Ionization Effects .....	139
Plasma Ionization Effects .....	140
Matching Overall Composition.....	141
Overcoming Physical Matrix Effects.....	142
What Is an Internal Standard?.....	143
Principles of Internal Standardization.....	144
Mass Effects and Internal Standardization .....	145
Typical Internal Standard Selections .....	146
Addition of Internal Standards.....	147
ISTD and Data Analysis Calculations .....	148
Selecting Internal Standards .....	149
Assigning an Internal Standard.....	151
Concentration of an Internal Standard .....	152
Reporting ISTD Results.....	153
What Is a Spectral Interference?.....	154
Elemental Isobaric Interferences.....	155

Molecular Polyatomic Spectral Interferences .....	156
Predictability of ICP-MS Interferences .....	157
Managing Spectral Interferences .....	158
How Large Is the Interference .....	159
Applying ORS Technology.....	160
Interference Correction Equations .....	161
Principle of Correction Equations.....	162
Correction Equation Dialog Boxes .....	164
Lead Equation Is a Unique Case.....	165
<b>OCTOPOLE REACTION SYSTEM (ORS).....</b>	<b>167</b>
In This Section .....	168
Octopole Reaction System Cell .....	169
Accessible Placement of the ORS.....	170
ORS Basic Principles.....	171
Helium: The Most Effective ORS Mode .....	172
Helium Cell Gas.....	173
Kinetic Energy Discrimination Effect .....	174
Principle behind Energy Discrimination.....	175
Bias Lens Setup for Efficient KED.....	176
Background Equivalent Concentration (BEC).....	177
Tuning He and H <sub>2</sub> ORS Gas Modes .....	178
Tuning Window for ORS Gas Tuning.....	179
Regular and High-Energy He Mode .....	180
Typical Settings: He and H <sub>2</sub> Modes.....	181
Analyte Sensitivity vs. He Flow Rate.....	182
Settings for He High-Energy Collision Mode .....	183
Hydrogen Reaction Gas Mode.....	184
Various Ion-H <sub>2</sub> Reactions and Rate Constants .....	185
Additional Ion-Molecule Reactions.....	186
Summary: Tuning of ORS He and H <sub>2</sub> Modes.....	187
Example: Improving Cr with He Mode .....	188
<b>MASSHUNTER ACQUISITION METHODS AND SEQUENCES FOR AGILENT 7700 ICP-MS SYSTEMS .....</b>	<b>189</b>
In This Section .....	190

Methods and Sequences .....	191
Necessity of a Method and Sequence File .....	192
Fundamentals of Acquisition .....	193
Peak Channels and Mass Sweeps .....	194
Executing Tune File Steps .....	195
Method Screens.....	196
Key Method Screens.....	197
Acquisition Modes .....	198
Acquisition Parameters: Spectrum Multi-Tune .....	200
Periodic Table and Mass Scale .....	201
Element Information .....	202
Acquisition Parameters: Tune and Time Settings.....	204
Peristaltic Pump Program for Autosampler .....	205
Advanced Peristaltic Pump Options .....	206
Saving and Running a Method.....	207
Acquisition Isotope Recommendations .....	208
Pre-Run Monitor for Manual Sampling.....	210
Creating a New Method: Method Wizard.....	211
MassHunter Sequences and Menu Selections.....	212
Edit Sample Log Table .....	214
Type: Keyword .....	224
Edit Sample Log Table .....	226
Sample Log Table Option 1: Simple Table Entry .....	227
Sample Log Table Option 2: Sequence Blocks .....	228
Example Blocks – Sequence Flow.....	230
Sample Log Table Option 3: Periodic Blocks .....	232
Whole List Display of Blocks.....	233
Running a Sequence.....	234
Chained Sequences .....	236
Chained Sequence Setup Screen.....	238
<b>DATA ANALYSIS METHODS AND PRINCIPLES OF QUANTIFICATION FOR FULLQUANT MODE .....</b>	<b>241</b>
In This Section .....	242
The Analytical Chemistry of IC-PMS .....	243

Data Analysis (DA) Module .....	244
Refining Raw ICP-MS Data .....	245
User Review Process.....	246
MassHunter DA Workflow .....	247
Create New Batch File .....	248
Load Batch File.....	249
Importing Files into DA Batch .....	250
Initial (Home) View of DA Batch .....	251
Toolbar Area of MassHunter DA .....	252
Creating Final ICP-MS Results .....	253
Editing Data Analysis: The DA Method Task Pane .....	254
Using Previously Configured Processing Details .....	255
Data Analysis Method.....	257
Analyte List Screen.....	258
FullQuant Window.....	259
FullQuant Outlier Screen .....	261
Report Templates and Worklist Actions.....	262
User Action: Process Batch Calculations .....	263
QC Check: Automatic Flagging of Outliers .....	265
Flags: Meanings in FullQuant Data Analysis .....	266
Batch Table Pane: Changing the Layout .....	267
On-Screen Reviewing of Results .....	268
Curve Pane Viewing .....	269
ICP-MS Results Reporting .....	270
Spectrum Pane for a Time Chart.....	271
Advanced ISTD Operations .....	273
Calibration.....	274
Viewing Calibration Curves .....	275
Common Calibration Curve Problems.....	276
Optimizing the Calibration Curve Fit .....	277
Calibration Curve Fitting Equation Options .....	278
Calibration Recommendations.....	280
Blanks and Solutions Used for Dilution .....	281

Application Figures of Merit .....	282
MassHunter Offline Data Analysis.....	283
<b>MASSHUNTER ANALYSIS CAPABILITIES .....</b>	<b>285</b>
In This Section .....	286
What Is SemiQuantitative Analysis? .....	287
SemiQuant Principles.....	289
SemiQuant: Acquisition Method Files.....	293
SemiQuant: Sequence Files .....	294
SemiQuant: Data Analysis Procedures .....	295
SemiQuant: Data Analysis Method Selection.....	296
SemiQuant: Initial View of Batch Results.....	297
SemiQuant: Response Factor Pane .....	298
Optimizing SemiQuant Results.....	299
SemiQuant: Updating the SQ Curve.....	301
Processing the SemiQuant Batch .....	302
SemiQuant: Reports .....	303
What Is Isotope Ratio Analysis?.....	304
Isotope Ratio Analysis: Acquisition Method.....	305
Isotope Ratios Data Analysis Setup.....	306
Isotope Ratio DA Method Setup.....	307
Isotope Ratio Batch Table.....	308
Isotope Ratio Reports.....	309
What Is Method of Standard Additions (MSA)? .....	310
Suitability of MSA: Pros and Cons.....	311
Editing Data Analysis for MSA .....	313
MSA Calibration and Result.....	314
MSA Results and Reports.....	315
<b>INTELLIGENT SEQUENCING FOR MASSHUNTER.....</b>	<b>317</b>
In This Section .....	318
Reviewing Run Performance .....	319
QA Major Divisions.....	320
Intelligent Sequencing Actions .....	322
QA Tuning Criteria .....	323

QC Database .....	328
QA Samples .....	330
Simulate Sequence .....	331
Data Analysis Method for Intelligent Sequencing.....	332
Processing Intelligent Sequence .....	333
View Intelligent Sequence Results .....	334
QC Database and ICP-MS Configuration.....	335
Main Database in Configuration.....	336
QC Main Parameters Setup.....	337
New QC Sample in Configuration .....	339
<b>AGILENT 7700 MAINTENANCE .....</b>	<b>341</b>
In This Section .....	342
Agilent ICP-MS Maintenance Video.....	343
Instrument Control Window .....	344
Maintenance LogBook.....	345
Maintenance Log .....	346
Running Time Feature .....	348
Maintenance Schedule .....	349
Sample Introduction Maintenance .....	351
Visual Aids for Maintenance .....	352
Torch Maintenance .....	353
Cone Maintenance .....	355
Extraction Lens Assembly Maintenance .....	356
Octopole Assembly Maintenance .....	357
Rotary Pump Maintenance.....	358
Multiplier Maintenance.....	359
Other Component Maintenance .....	360
Hydrogen Generator.....	362
<b>INTRODUCTION TO HYPHENATED ICP-MS TECHNIQUES.....</b>	<b>363</b>
In This Section .....	364
Features of Hyphenated Techniques.....	365
Laser Ablation-ICP-MS.....	366
Speciation.....	374

Complementary Techniques .....	376
ICP-MS Interfacing Options .....	377
GC-ICP-MS .....	378
HPLC-ICP-MS .....	381
Capillary Electrophoresis-ICP-MS .....	384
Summary .....	386
<b>ICP-MS RESOURCES .....</b>	<b>387</b>
In This Section .....	388
Staying Up-to-Date .....	389
ICP-MS Website .....	390
ICP-MS Information CD .....	391
The Agilent ICP-MS Journal .....	392
E-Seminars .....	393

# Agilent 7700 Series ICP-MS Techniques and Operation

Course Number R1777A  
Student Laboratory Manual  
Appendix





**Agilent Technologies**

# Agilent 7700 Series ICP-MS Techniques and Operation

Course Number R1777A

Student Laboratory Manual and  
Appendix

Agilent 7700 MassHunter Workstation  
Printed in September, 2009

## **Notice**

The information contained in this document is subject to change without notice.

Agilent Technologies makes no warranty of any kind with regard to this material, including but not limited to the implied warranties of merchantability and fitness for a particular purpose.

Agilent Technologies shall not be liable for errors contained herein or for incidental, or consequential damages in connection with the furnishing, performance, or use of this material.

No part of this document may be photocopied or reproduced, or translated to another program language without the prior written consent of Agilent Technologies, Inc.

Agilent Technologies, Inc  
3750 Brookside Pkwy  
Suite 100  
Alpharetta, GA 30022

© 2009 by Agilent Technologies, Inc.

All rights reserved

Printed in the United States of America

## Table of Contents

<b>LABORATORY EXERCISE: AGILENT 7700 ICP-MS START UP AND TUNING .....</b>	<b>1</b>
In this laboratory exercise you will:.....	2
Requirements .....	3
Instrument Start Up.....	4
Tuning: Normal mode.....	5
Tuning: He Mode.....	7
Tuning: High Energy He Mode .....	8
Tuning: H <sub>2</sub> Mode.....	9
<b>LABORATORY EXERCISE: ACQUIRE A SIMPLE MASS SPECTRUM.....</b>	<b>11</b>
Requirements .....	13
Acquire Data.....	14
Review Data.....	16
<b>LABORATORY EXERCISE: MULTI-TUNE METHOD AND SEQUENCE.....</b>	<b>19</b>
Requirements .....	21
Method Parameters .....	22
Sequence Set up and Run.....	26
Method and Sequence Review and Conclusions .....	28
<b>LABORATORY EXERCISE: DATA ANALYSIS SET UP AND EVALUATION .....</b>	<b>29</b>
Requirements .....	31
Create Batch, Import Data and build DA Method .....	32
Evaluate results .....	34
Outliers.....	35
Report.....	36
Data Analysis Conclusions .....	37
<b>APPENDIX: TIME RESOLVED ANALYSIS, TRANSIENT SIGNALS AND MASSHUNTER SOFTWARE .....</b>	<b>39</b>
In This Section .....	40
Time Resolved Analysis .....	41
TRA Parameters.....	42
TRA Parameters.....	43
TRA Parameters.....	44
MassHunter Chromatographic Data Analysis .....	46

MassHunter Chromatographic Data Analysis .....	47
MassHunter Chromatographic Data Analysis .....	48
Compound Independent Calibration .....	49
Compound Independent Calibration .....	50
<b>APPENDIX: ANALYSIS OF TRACE METALS IN ORGANIC SAMPLES .....</b>	<b>51</b>
In This Section .....	52
Organic Samples .....	53
Organic Samples .....	54
Practical Considerations.....	55
Practical Considerations.....	56
Organic Solvent Configuration .....	57
Hardware Recommendations .....	58
Hardware Recommendations .....	59
Oxygen Addition.....	60
Oxygen Addition.....	61
Oxygen Addition.....	62
Spectroscopic Challenges .....	63
Method of Standard Additions.....	64
Summary .....	65
User Forum Resources.....	66
<b>APPENDIX: INTEGRATED SAMPLE INTRODUCTION SYSTEM.....</b>	<b>67</b>
In This Section .....	68
Integrated Sample Introduction System.....	69
ISIS Applications .....	70
Hardware Configurations.....	71
Hardware Configurations.....	72
Constant Flow Nebulization .....	73
Constant Flow Nebulization .....	74
On-line Auto Dilution.....	75
On-Line Auto Dilution.....	76
On-line Auto Dilution.....	77
Discrete Sampling.....	78
Discrete Sampling.....	79

Discrete Sampling.....	80
Hydride Generation.....	81
Hydride Generation.....	82
<b>APPENDIX: ULTRA-TRACE LEVEL ULTRA-HIGH PURITY APPLICATIONS .....</b>	<b>83</b>
In This Section .....	84
Common Sources of Contamination.....	85
Sample Preparation Environment .....	86
Laminar Flow Hood.....	87
Laminar Flow Hood.....	88
The Operator .....	89
Deionized Water .....	90
Labware.....	91
Labware.....	92
Labware.....	93
Pipette Tips .....	94
Cleaning Tanks .....	95
Cleaning Procedures .....	96
Cleaning Procedures .....	97
Cleaning Procedures .....	98
Acid Redistillation .....	99
Concentrated Acids and Reagents .....	100
Certified High Purity Chemicals.....	101
Preventing Chemical Contamination .....	102
Preventing Chemical Contamination .....	103
Sample Preparation .....	104
Useful Laboratory Items .....	105
Conclusions.....	106